




# Fitting Instructions

Disposable respirators are only most effective when there is a good seal between the edges of the respirator and your face. The instant this seal is broken, protection is compromised as contaminated air can leak in through any gaps.

**These fitting instructions must be followed each time a 3M™ Aura™ 9400+ Series respirator is worn.**



Make sure that your face is clean shaven. Respirators should not be worn with stubble, beards or other facial hair under the area of the face seal as these can prevent a good seal to the face. Make sure that long hair is tied back and jewellery is removed so that it does not interfere with the seal to the face.

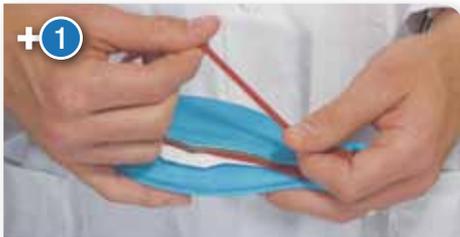

**1.** Pre-stretch around entire length of each strap by pulling at 3cm intervals between both hands.



**2.** With the reverse side up and using the tab, separate the top and bottom panels of the respirator to form a cup shape. Bend slightly at the centre of the noseclip.

**3.** Ensure that both panels are fully unfolded.



**4a.** Cup the respirator in one hand with the open side towards your face.



**4b.** Take both straps in your other hand. Hold the respirator under your chin, with the nosepiece facing upwards and pull the straps over your head.



**5.** Locate the upper strap across the crown of the head and the lower strap below your ears. The straps must not be twisted. Adjust the top and bottom panels for a comfortable fit, ensuring that the panels and tab are not folded in.



**6.** Using both hands, mould the noseclip to the shape of the nose to ensure a close fit and a good seal.

**▲ The respirator may not fit as well if you pinch the noseclip using one hand. Use two hands.**



**7.** Perform a fit-check by covering the front of the respirator with both hands taking care not to disturb its fit. **Inhale** sharply.

If air leaks around the nose, readjust the noseclip to eliminate leakage then repeat the fit check. If air leaks at the respirator edges, work the straps back along the sides of the head to eliminate leakage then repeat the fit check. **If you cannot achieve a proper fit, DO NOT enter the hazardous area. Consult your supervisor.**

The Power to Protect Your World™



**Safety Solutions  
3M United Kingdom plc**  
3M Centre  
Cain Road, Bracknell  
Berkshire RG12 8HT  
Tel: 0870 60 800 60  
www.3M.co.uk/ohes

**Safety Solutions 3M  
Ireland Limited**  
The Iveagh Building  
The Park  
Carrickmines  
Dublin 18  
Tel: 1 800 320 500



Distributed by Industrial Scientific Supplies, a division of Camlab Ltd

Camlab Ltd, Norman Way Industrial Estate, Over, Cambridge, CB24 5WE  
Telephone: +44 (0)1954 233145  
Email: sales@camlab.co.uk  
www.industrialscientific.co.uk



Scan this QR code with your smartphone to be directed to further information on 3M™ Aura™ 9400+ Disposable Respirators

14048  
Please recycle. Printed in UK.  
© 3M 2011. All rights reserved.